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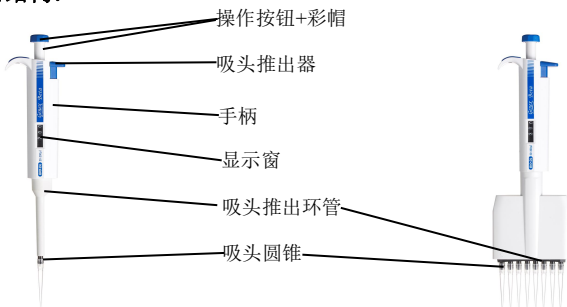
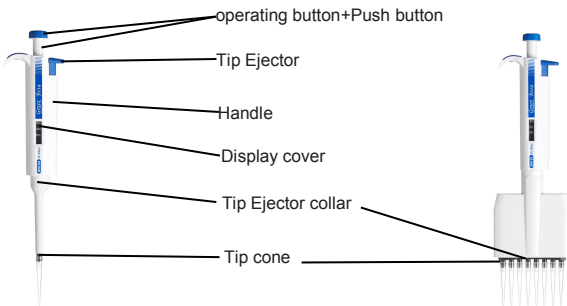
2、Pipette Operation:

A、Volume setting

*The volume of the pipettor is clearly shown through the handle grip window. The delivery volume (variable volume pipettors only) is set by turning the thumb button clockwise or anticlockwise .

*When setting the volume. The desired delivery volume clicks into place. The digits are completely visible in the display window. The selected volume is within the pipettor's specified range.

B、Pipetting technique

Forward pipetting:

*Make sure that the tip is firmly attached to the tip cone. For best results the thumb button should be operated slowly and smoothly at all times, particularly with viscous liquids.

*Hold the pipettor vertically during aspiration. Make sure that the liquid and container vessel are clean and that the pipettor, tips and the liquid are at the same temperature.

1) Depress the operating button to the first stop.

2) Place the tip(s) just under the surface of the liquid (2-3 mm) and smoothly release the operating button. Carefully withdraw the tip from the liquid, touching against the edge of the container to remove excess.

3) Liquid is dispensed by gently depressing the thumb button to the first stop. After a short delay continue to depress the thumb button to the second stop. This procedure will empty the tip(s) and ensure accurate delivery.

4) Release the thumb button to the ready position. If necessary change the tip(s) and continue with pipetting.

3、STORAGE:

*When not in use it is recommended that your pipettor is stored in a vertical position.

Shanghai BIO-DL Science Instrument Co., Ltd.
Building 12, No.201 of Minyi road, Songjiang district,
Shanghai, China 201101.

Tel: 021—37012760

Fax: 021—37012768

<http://www.bio-dl.com.cn>

Please scan the QR code to get more instructions